

Stationary Basket Installation and Setup

- 1) Set the Paddle Height to 25 mm \pm 2 mm from the bottom of the vessel. [SEE FIGURE 1](#)
 - 2) Assemble the holder loosely to the cover with nuts all the way at the bottom and the top of the thread. [SEE FIGURE 2](#)
 - 3) Install the 1 cm spacer (DEPSET-SB) on the guide rod. If a 25 mm spacer is already installed, install the 1 cm spacer on top. [SEE FIGURE 3](#)
 - 4) Install the cover assembly in the vessel and rotate the paddle so that it is directly beneath the basket and touching the basket. [SEE FIGURE 4](#)
 - 5) Lower the top nuts so they touch the cover. [SEE FIGURE 4](#)
 - 6) Raise the drive head, remove the cover assembly and tighten bottom nuts to cover. The height is now set. [SEE FIGURE 5](#)
 - 7) Remove spacer. [SEE FIGURE 6](#)
- For DEPSET-SB continue to step 3.
For DEPSET SBA see below.
- a) Install the DEPSET-SBA spacer as per [FIGURE 3a](#).
 - b) Install the cover assembly in the vessel so the basket touches the DEPSET-SBA. [SEE FIGURE 3b](#)
 - c) Lower the top nut so it touches the cover. [SEE FIGURE 4](#)
 - d) Lift the Drive head and remove the spacer. Remove the cover and tighten the bottom nuts to cover. [SEE FIGURE 5](#)
The height is now set.



FIGURE 1

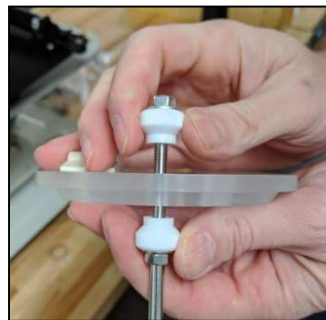


FIGURE 2



FIGURE 3



FIGURE 3a



FIGURE 3b

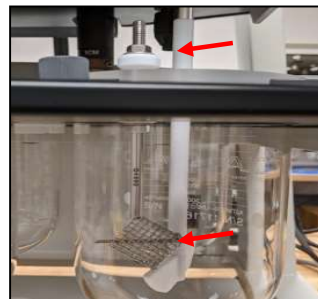


FIGURE 4

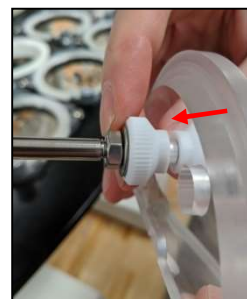


FIGURE 5



FIGURE 6